


PROTOCOL

BetterBuffer
Sequence Enhance/Dilution Buffer

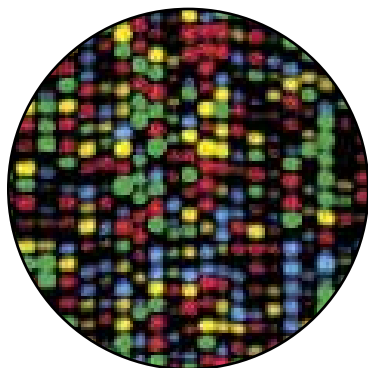

Part Number: DAF-XX
Version: 4.0
Storage: Store @ -20°C
Related Accessories: microCLEAN, MAP100C 96 Well AMP plates,
10X CE Running Buffers

- Dilute the ABI BigDye Mix with BetterBuffer to achieve the desired concentration*
- Mix by pipetting
- Sequence according to protocol
- Remove excess dye

*1 in 6, 1 in 12, 1 in 15 or higher. The choice of dilution is dependent on size, quantity and quality of the DNA template.

BigDye is a registered trademark of Applied Biosystems

Supplied as a clear solution in multiple of 0.3, 1.5, 25 & 250 ml aliquots
For Research Only